

Software installation

1. Installation of system program

- (1) Turn on the power switch of a personal computer and start Windows[®].
- (2) Exit other programs.
- (3) Set a system program CD-ROM to the drive. The system setup window is displayed.
- (4) Install system program by the window indication.
- (5) After the setup, take out the system program CD-ROM.

2. Installation of Microsoft[®] Excel

This software is required when displaying measurement data at worksheet by using Excel Macro.

- (1) Turn on the power switch of a personal computer and start Windows[®].
- (2) Exit other programs.
- (3) Set a Microsoft Excel CD-ROM to the drive. The system setup window is displayed.
- (4) Operate Microsoft[®] Excel by the window indication. Select "Standard" for installation.
- (5) After the setup, take out the Microsoft[®] Excel CD-ROM and restart the personal a computer.

* For Excel 2007, change the macro setting by the following procedures.
(Without this setting, the error message is displayed and macro can't be started.)

1. Start Microsoft[®] Excel. Click [Office] button of the upper left and [Excel Option (I)] button of the bottom.
2. Click "Security center" of the left and [Security center setting (T)] button in "Excel Option" window.
3. Click "Macro setting" of the left, "Enable all macros (E)", and [OK] button in "Security center" window.
4. Click [OK] button in "Excel Option" window.

3. Installation of printer drivers

Previously install printer drivers into PCs to install NSX-100 PC software.

If you run print or preview without installing printer drivers, "Run-time error '380'" occurs.

4. Uninstallation

When the program does not operate properly, uninstall the system program and install it again.

- (1) Click [Start] button, "Setup" and "Control Panel".
- (2) Click "Application Addition and Deletion."
- (3) Select a program to delete from the list of "Installation and Deletion".
- (4) Click "Addition and Deletion" button. Uninstallation window is displayed.
- (5) Select "Automatic Uninstallation" and click [Next] button. Uninstall the system program.